

SWAB SYSTEMS for environmental monitoring in the Pharmaceutical Industry

Ready and easy to use, Copan swabs are now available from Oxoid. These individually sterile-wrapped swabs which are world-renowned for quality and excellence, have been developed by experts in the pre-analytical phase of microbiology:

Collection and transportation of microbiological samples

- Ideal for use in the monitoring of clean rooms, isolators and other critical areas.

Single tube systems

- For sampling and immediate streaking onto appropriate culture medium.

Sterility Assured

- Irradiated at 25K Gy for sterility assurance.
- Vaporized Hydrogen Peroxide (VHP) resistant packaging.
- Triple wrapped for use in critical areas such as clean rooms and isolators.

Easy to Use and Store

- No glass or sharp edges.
- No refrigeration required, store at room temperature (5 - 25°C).
- Long shelf-life for prolonged storage.

Effective

- Inactivates common cleaning agents and dissolves bio-films.

Quality Certification

- Quality certificates and technical files available on request to satisfy quality assurance requirements.







Procedure

Single tube kit - for immediate streaking onto culture plate

1. Peel open pack.
2. Remove tube cap.

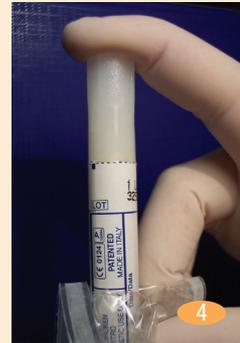

3. Insert swab in tube.
4. Close cap.
5. Squeeze base of tube to moisten swab tip.

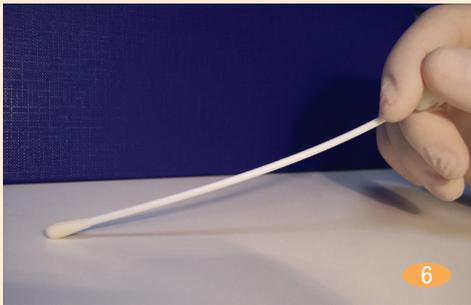

6. Sample surface or equipment.
7. Roll swab on appropriate culture medium.

Ordering information

Product Description	Pack Size	Order Code
Triple Wrapped	200	TS0954C



DEDICATED TO MICROBIOLOGY

www.oxid.com

Oxoid Ltd, Wade Road, Basingstoke, Hants, RG24 8PW, UK. Tel: +44 (0) 1256 841144 Fax: +44 (0) 1256 463388 Email: oxid@oxid.com

